

IMAGING AND ANALYSIS SOFTWARE


Gel/Blot Alignment Templates for the ChemiDoc™ and ChemiDoc MP Imaging Systems

Quick Guide


The Bio-Rad gel/blot alignment templates enable you to quickly and consistently place gels and blots on the imaging systems' sample trays.

The Gel/Blot Alignment Templates (catalog #12003915) contain the following:

- Alignment template for small gels, 9.5 x 7.6 cm

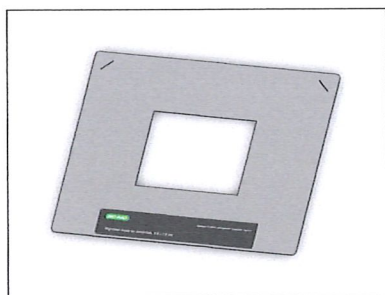

- Alignment template for medium gels, 15.5 x 12.4 cm

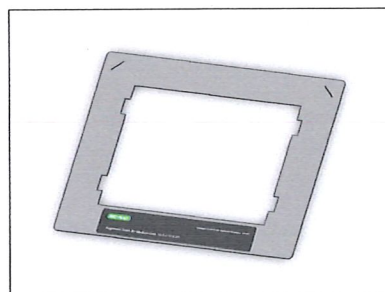

You can use the gel/blot alignment templates with all supported tray types. For more information about supported tray types, see the ChemiDoc and ChemiDoc MP Imaging Systems with Image Lab™ Touch Software User Guide (#10000062126).

Prepare the Gel/Blot Alignment Template

- Select the gel/blot alignment template that matches the size of your sample.
- Bend a corner at the pre-cut slit to make a tab. Use the tab to remove the template from the tray after use.
- Place the gel/blot alignment template on the sample tray.
- Place the gel/blot or gel/blot tray on the open area of the template.



Alignment template for medium gels

Clean and Store Templates

It is important to keep the templates clean and to store them flat.

- After using a gel/blot alignment template, rinse it with water and allow it to air dry.
- Store the gel/blot alignment templates flat to prevent prevent warping.

BIO-RAD



**Bio-Rad
Laboratories, Inc.**

Life Science
Group

Web site bio-rad.com USA 1 800 424 6723 Australia 61 2 9914 2800 Austria 43 1 877 89 01 177 Belgium 32 (0)3 710 53 00 Brazil 55 11 3065 7550
Canada 1 905 364 3435 China 86 21 6169 8500 Czech Republic 420 241 430 532 Denmark 45 44 52 10 00 Finland 358 09 804 22 00
France 33 01 47 95 69 65 Germany 49 89 31 884 0 Hong Kong 852 2789 3300 Hungary 36 1 459 6100 India 91 124 4029300
Israel 972 03 963 6050 Italy 39 02 216091 Japan 81 3 6361 7000 Korea 82 2 3473 4460 Mexico 52 555 488 7670 The Netherlands 31 (0)318 540 666
New Zealand 64 9 415 2280 Norway 47 23 38 41 30 Poland 48 22 331 99 99 Portugal 351 21 472 7700 Russia 7 495 721 14 04
Singapore 65 6415 3188 South Africa 27 (0) 861 246 723 Spain 34 91 590 5200 Sweden 46 08 555 12700 Switzerland 41 026674 55 05
Taiwan 886 2 2578 7189 Thailand 66 662 651 8311 United Arab Emirates 971 4 8187300 United Kingdom 44 020 8328 2000

